

Qiagen Gel Extraction

Rationale:	
Special Observations:	
Results:	
Interpretation:	

Experiment Date:	Source: Caleb Radens,
Experiment Time	Ali Awan
Primary Experimenter (contact):	Assembled: 7/19/2012
Other Experimenters:	

DNA details	~ug DNA in gel	Yield (ng/uL)

Procedure:

Critical Steps:

- Minimize UV exposure to DNA and yourself
 - Protect your eyes from the UV light
- Turn on 50 dc water bath**
- Stain gel with ethidium bromide**
- Approximately 1 hour, or just enough to visualize ladder and DNA of interest
- Visualize gel to check DNA band of interest is present and take a picture (See: Weill Gel Pictures)**
- Be sure to bring or memorize the ladder bands and bring documentation of what DNA is in which lane
 - Minimize UV exposure time
- After taking picture, use a razor blade to cut the DNA fragment of interest out of the gel**

- Minimize UV exposure time
 - Cut as close to the DNA as possible
 - After making four incisions around band(s) turn off the UV then excise your fragment.
Flip the fragment on its side, turn on UV, then cut off any extra non-glowing gel
 - You want to have a slice weighing near of less than 200 mg
- Put the fragment into a pre-weighed and labeled 2 mL centrifuge tube**
 - Determine the weight of the fragment**
 - Add (3)x(Gel fragment weight in mg) uL of buffer QG to tube**
 - Ex) 0.202 g gel slice → add 606 uL buffer QG
 - Vortex tube, then incubate in 50 dc water bath for at least 20 minutes, vortexing every couple of minutes**
 - At this time, label Qiaquick spin column(s)
 - Transfer dissolved gel sample to the top of a labeled Qiaquick spin column**
 - Centrifuge @ 13,000 rpm 1 minute, discard flow-through**
 - Add 500 uL of buffer QG to column**
 - Centrifuge column @ 13,000 rpm 1 minute, discard flow-through**
 - Add 750 uL buffer PE to column**
 - Start pre-warming buffer EB to 70 dc**
 - Centrifuge column @ 13,000 rpm 1 minute, discard flow-through**
 - Add 750 uL buffer PE to column (Yes, again)**
 - Centrifuge column @ 13,000 rpm 1 minute, discard flow-through**
 - Be sure to vigorously shake out the residual liquid in the flow through tube
 - Centrifuge column with the lid open for 2 minutes @ 13,000 rpm**
 - Transfer spin column to labeled tube**
 - Add 30 uL of buffer EB directly to center of membrane**
 - It's ok to touch the membrane to dispel a residual droplet is on pipette tip
 - Wait AT LEAST 5 minutes (with lid closed)**
 - May wait longer if you're patient, but don't wait more than 30 or so minutes
 - Centrifuge column @ 13,000 rpm 1 minute, if there is any leftover EB visibly on the membrane, rotate the column 180 degrees and centrifuge again**
 - Expect very low concentration of DNA (~10 – 30 ng/uL from 1 ug DNA)**